rPLD (CLA Antigen)

Vial of 100µg / Flacon de 100µg

Ref. VP003A

STORE AT
2-8°C

For research use only

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CONSERVER A
2-8°C

Origin:

rPLD (CLA Antigen), highly purified and stabilised. rPLD is a recombinant form of phospholipase D, an important, conserved *C. pseudotuberculosis* bacterial virulence factor. rPLD was identified as an antigen for IgG antibodies developed in the caseous lymphadenitis (CLA) disease in sheep or goat sera.

Presentation:

Vial containing approximately 100 µg of rPLD lyophilised in presence of additives and preservatives.

Reconstitution:

Each vial must be restored with 0.5 ml distilled water in order to obtain a preparation at 200µg/ ml. It can then be diluted using the appropriate solution.

Excipients:

– Phosphate, Mannitol.

Purity:

Highly purified recombinant PLD, with a major band of about 31 KDaltons on SDS-PAGE. Presence of Dimers can occur.

Safety

Any product of biological origin must be handled with all the required cautions, as being potentially infectious.

Use:

For *in vitro* use only.
All research studies and protocols where a source of highly purified rPLD is necessary.

Stability:

Restored :

- 24 hours at room temperature
- 72 hours at 2-8°C
- 2 months at -20°C or below

Lyophilized : - Until the expiration date printed on the vial.

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CONSERVER A
2-8 °C

ANALYSIS CERTIFICATE

ANALYTICAL DATA	SPECIFICATIONS
1. Protein Content µg/vial	> 90 µg/vial
2. SDS-PAGE (12 % acrylamide) 1 major band of about daltons	1 major band of about 31,000 daltons

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DATE :

CONCLUSIONS : Passed/Conforme Refused/Non-conforme

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